Chloramphenicol resistant pUC Vector

pHSG398 DNA

Code No. 3398

Size : 25 μ g (0.5 OD)

Shipping at – 20°C

Stored at – 20°C

* 2 years from date of receipt under proper storage conditions.

Lot No.

Concentration :

μ g/ μ l

Volume :

μ l

Description:

pHSG398 is a pUC type of cloning vector plasmid possessing chloramphenicol resistancy. This plasmid is suitable for DNA sequencing by dideoxy method. It may allow successful cloning of DNA fragment longer than the case of M13 phage vector used. Since it has multi-cloning sites on *lacZ* region, any DNA insertion into this plasmid can be easily verified using plates containing IPTG and X-Gal. Moreover, expression of the cloned DNA can be performed by using of *lac* promoter. M13 primers are available for DNA sequencing.

Using Deletion Kit for Kilo-Sequencing (Cat. #6030), DNA sequencing of the cloned DNA with kilo-bases can also be performed.


pHSG398 contains the consensus multi-cloning sites as same as pUC18.

Form: 1 mM EDTA
10 mM Tris-HCl, pH 8.0

Preparation: cccDNA was purified by CsCl-EtBr ultracentrifugation

Base pairs: 2,227 bp

Multicloning site of pHSG398:

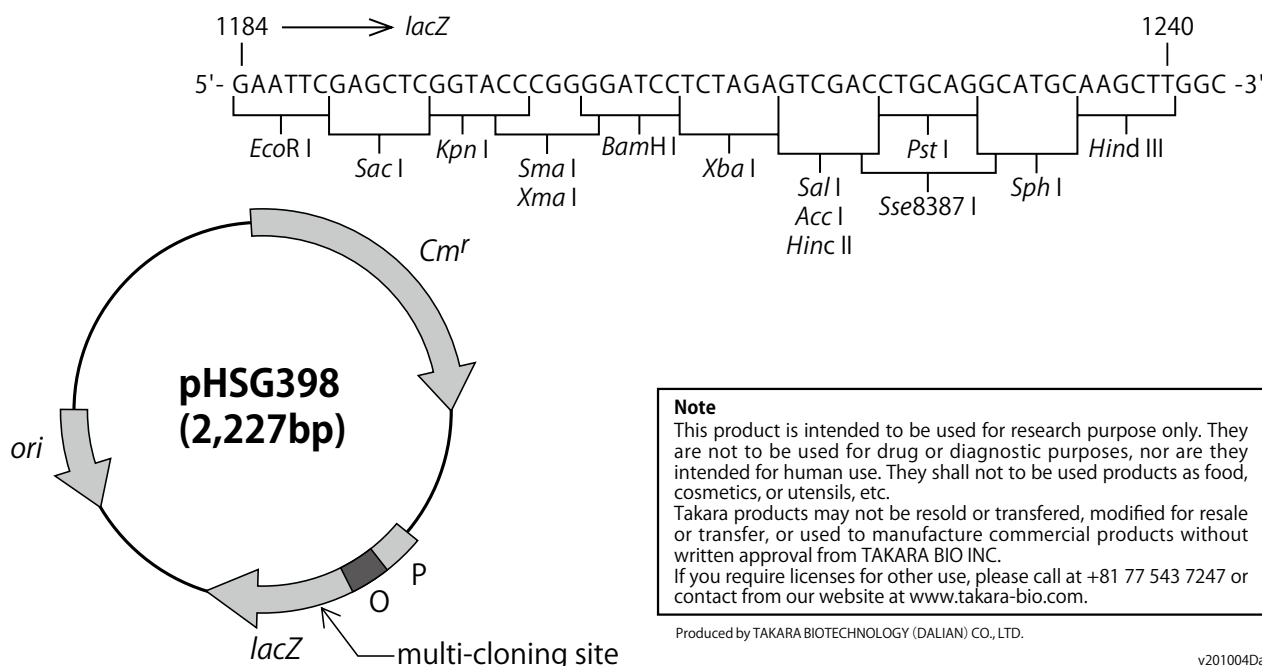Purity:

1. Contains over 70% double-stranded covalently closed circular form I (RF I) DNA.
2. Shown to retain cloning site by dideoxy sequencing method.
3. Shown to be cleaved at a single site by restriction enzymes *EcoR* I, *Sac* I, *Kpn* I, *Sma* I, *Bam* H I, *Xba* I, *Sal* I, *Pst* I, *Sph* I and *Hind* III.

Usage:

- Cloning of a target gene and its expression using *lac* promoter
- DNA sequencing by using M13 primers
- Determination of long DNA sequences using Deletion Kit for Kilo-Sequencing (Cat. #6030)

Reference:

Takeshita, S., Sato, M., Toba, M., Masahashi, W. and Hashimoto-Gotoh, T. (1987) *Gene* **61**, 63-74.

Note

This product is intended to be used for research purpose only. They are not to be used for drug or diagnostic purposes, nor are they intended for human use. They shall not to be used products as food, cosmetics, or utensils, etc.

Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from TAKARA BIO INC.

If you require licenses for other use, please call at +81 77 543 7247 or contact from our website at www.takara-bio.com.

Produced by TAKARA BIOTECHNOLOGY (DALIAN) CO., LTD.

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