## RNase A

### 1 Contents

RNase A		
REF	740505	740505.50
RNase A (lyophilized)	100 mg	50 mg

### 2 Product description

The RNase A is primarily intended to be used with NucleoSpin® and NucleoBond® kits. See individual protocols for detailed information.

However, if the RNase A is used for another than the recommended purpose performance may differ considerably from its expected one.

### 3 RNase A specifications

- Source:** Bovine pancreas  
**Form:** 100 mg / or 50 mg white freeze-dried material  
**DNase activity:** Not detectable if used in above mentioned kits

### 4 Storage conditions and preparation of working solution

Lyophilized RNase A can be stored at room temperature (18–25 °C) for at least one year.

A: Dissolved in water: Store the RNase A solution at 4 °C for up to 3 months. For longer storage, the RNase A solution should be divided into small aliquots and stored at -20 °C.

B: Dissolved in Resuspension Buffer A1/S1/S1-EF/RES/RES-EF/RES-BAC: Buffer containing RNase A can be stored at 4 °C for at least 6 months.

RNase A		
REF	740505	740505.50
RNase A (lyophilized)	100 mg	50 mg
	Dissolve in H <sub>2</sub> O or directly in Resuspension Buffer A1 / S1 / S1-EF / RES / RES-EF / RES-BAC*	

\* Please refer to the individual protocols for detailed information.

## 5 Safety instructions

RNase A contains hazardous contents. Wear gloves and goggles!

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### 5.2 GHS classification

Only harmful features must not be labeled with H and P phrases until 125 mL or 125 g.

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