pT7-IRES His-C DNA

Code No. 3291

Size: 20 µg

Conc.: 0.5 µg/µl

Chain length: 3,429 bp

Quality Control Data:


Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

* 2 years from date of receipt under proper storage conditions.

Usage:

Protein expression using Human Cell-Free Protein Expression System

Vector map for pT7-IRES His-C DNA:

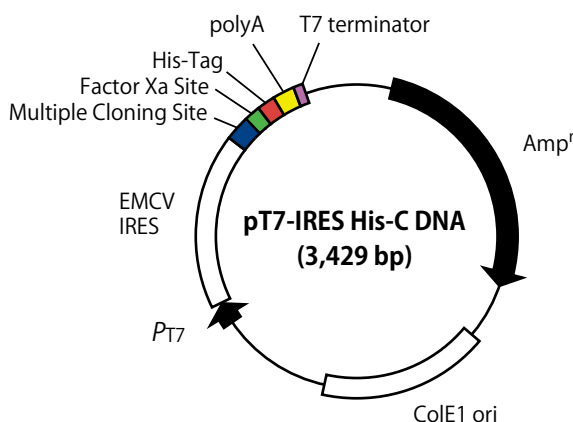Regarding protocol for protein synthesis, please refer to the product manual for Human Cell-Free Protein Expression System (Cat. #3281).

Description:

pT7-IRES DNA series are expression vectors designed for Human Cell-Free Protein Expression System. Tag sequence such as His-Tag or c-Myc Tag, Factor Xa cleavage site, Multiple Cloning Site (MCS), polyA, and T7 terminator are located at the downstream of T7 promoter and EMCV IRES. There are vectors with some arrangements for the kind of Tag and the location.

pT7-IRES His-C DNA is an expression vector including His-Tag sequence. Using pT7-IRES His-C DNA together with Human Cell-Free Protein Expression System (Cat. #3281) enables the synthesis of your target protein as C-terminal fusion of His-Tag. Factor Xa cleavage site is inserted so that His-Tag can be removed from the synthesized fusion protein.

Target gene cloned into MCS in frame is transcribed as RNA-containing-EMCV IRES under the control of T7 promoter. By the effect of EMCV IRES, designed to promote protein translation initiation, efficient high level protein synthesis can be performed in Human Cell-Free Protein Expression System.

Form: 10 mM Tris-HCl, pH 8.0
1 mM EDTA

Storage: -20°C

Preparation: Purified by ion-exchange column.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at www.takarabio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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		<u>Nco I</u>																	
<u>EMCV IRES</u>																			
5'-TAACGT.....TAATATGGCCACAACC	ATG	GCT	AGC	CAT	ATG	GAG	CTC	CTC	GAG	GGA	TCC	GAA	TTC	ACT	AGT	GTC	GAC	CTG	CAG
3'-ATTGCA.....ATTATACCGGTGTTGG	TAC	CGA	TCG	GTA	TAC	CTC	GAG	GAG	CTC	CCT	AGG	CTT	AAG	TGA	TCA	CAG	CTG	GAC	GTC
		Met																	
		Ala	Ser	His	Met	Glu	Leu	Leu	Glu	Gly	Ser	Glu	Phe	Thr	Ser	Val	Asp	Leu	Gln
		<u>Xba I</u>																	
<u>Factor Xa</u>																			
		<u>His-Tag</u>																	
		<u>End</u>																	
TCT	AGA	ATC	GAA	GGG	CGC	CAC	CAT	CAC	CAT	CAC	CAT	TAG	GTAATC-3'						
AGA	TCT	TAG	CTT	CCC	GCG	GTG	GTA	GTG	GTA	GTG	GTA	ATC	CATTAG-5'						
Ser	Arg	Ile	Glu	Gly	Arg	His	His	His	His	His	His	His							

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