CLEARLY Protein Ladder (Unstained)

Code No. 3453A Size: 500 μ l

Description :

CLEARLY Protein Ladder (Unstained) is a mixture of 11 recombinant highly purified proteins. When this ladder is analyzed by SDS-PAGE and stained with Coomassie Brilliant Blue R-250, it is resolved into quite clearly identifiable bands from 10 to 250 kDa. This ladder is intended for use as a precise size standard when performing SDS-PAGE to calculate the molecular weight of a protein of interest. The 30 and 80 kDa bands have quadruple the mass of the other bands and serve as reference indicators. This ladder is pre-mixed with loading buffer and ready to load. Pre-heating is not required. 5 μ l of this ladder is loaded per lane of SDS-PAGE minigel for standard use. In this case, this ladder is sufficient for 100 lanes.

Storage: -20°C

- * This ladder is stable for 6 months if stored at 4°C.
- * This ladder is stable after 50 times of freeze-thaw cycles.


Application Example :

1. Thaw CLEARLY Protein Ladder (Unstained) at room temperature. Confirm the ladder is completely dissolved. Pre-heating is not required.
2. Load 5 μ l per lane of SDS-PAGE minigel. Run SDS-polyacrylamide gel electrophoresis.
3. Perform staining with Coomassie Brilliant Blue R-250, etc.

Note :

It is recommended to use 5 - 20% gradient SDS-polyacrylamide gel for electrophoresis. If the concentration of gel is too low, the mobility of bands of low molecular weight is faster than the mobility of BPB. If the concentration of gel is too high, high molecular weight bands may not separate well or they may not migrate.

< Electrophoresis Result (12% SDS-PAGE) >

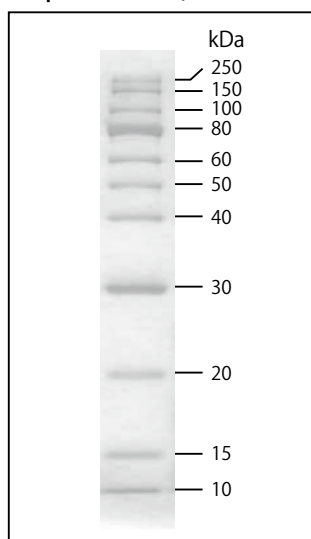Staining with LabSafe GEL Blue (Cat. #786-35)

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc. If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com. Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements. All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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