β -globin (human) Primer set

Code No. 3868

Size: 500 pmol each

Description :

This product is a set of PCR primers for amplification of β -globin region of human genome DNA. By combining two of these primers, nine different lengths of DNA fragment can be amplified. Therefore, these primers can be used as a positive control in PCR reaction when confirming whether template DNA is appropriate as PCR template.

Components : each 500 pmol

1. PC03 (forward) d (ACACA ACTGT GTTCA CTAGC)
2. PC04 (reverse) d (CAACT TCATC CACGT TCACC)
3. GH20 (forward) d (GAAGA GCCAA GGACA GGTAC)
4. GH21 (reverse) d (GGAAA ATAGA CCAAT AGGCA G)
5. KM29 (forward) d (GGTTG GCCAA TCTAC TCCCA GG)
6. KM38 (reverse) d (TGGTC TCCTT AAACC TGTCT TG)

Form : Lypophilized

Storage : -20°C

Production Method : Produced by Solid-phase synthesis method.

Quality Control Data :

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Reconstitution:

Reconstitute by adding sterile purified water or TE Buffer (pH 7.5 - 8.0). Reconstituted solution should be kept at -20°C.

Amplified fragment (bp) :

		reverse		
		PC04	GH21	KM38
forward	PC03	110	250	167
	GH20	268	408	325
	KM29	205	345	262

PCR condition :

94°C	1 min	} 25 - 35 cycles
55°C	2 min	
72°C	1 min	

Reference :

Saiki R K, Gelfand D H, Stoffel S, Scharf S J, Higuchi R, Horn G T, Mullis K B, and Erlich H A. *Science*. (1988) **239**: 487-491.

Note :

Primer PC03 has the same sequence as described in the reference, however, some template DNA might not be amplified efficiently using this primer. In this case, it is advisable to use other primer pair.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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