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The screenshot shows the Exiqon website header with the logo and navigation links: Home, Search, Contact, Print, Basket, and Sign In. The Sign In link is circled in red and labeled with a circled '1'. Below the header is a navigation bar with links: Products, Services, Applications, Resource Center, Ordering, About Exiqon, and Investor. The main content area is titled 'Sign in' and contains a form with the following elements: a 'Username or email:' input field, a 'Password:' input field, a 'Remember me next time.' checkbox, a 'Sign in' button, and two links: 'Register' (circled in red and labeled with a circled '2') and 'Forgot password?'.

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ExiLERA™ qPCR primers

This web tool is used for the design of custom LNA™-enhanced qPCR primer sets for mRNA and lncRNA from any organism. The algorithm is optimized for design of primers for RNA sequences of more than 55 nucleotides. For shorter sequences please refer to our [microRNA qPCR design tool](#)

Related products

The LNA™ primer sets designed using this software are optimized for use with the following products:

→ [cDNA Synthesis and SYBR® Green Master Mix Kits](#)

The screenshot shows the design interface with the following elements circled in red and numbered:

- ①: Text input field for "Enter a name for your project".
- ②: Dropdown menu for "Select organism" with "Human" selected.
- ③: Text input field for "Specify your target(s)" with the placeholder "By sequence ID (GenBank, RefSeq, Ensembl or keyword e.g. NM_001101)".
- ④: Large text area for "By Sequence (fasta, genbank or embi)".
- ⑤: "Next" button.
- ⑥: "Advanced options" button.

Other visible elements include a "Clear data" button, a "Send" button, and a checkbox for "Automatically select intron spanning assays".

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The screenshot shows the progress screen with the following elements:

- EXIQON logo with tagline "Seek Find Verify".
- Progress indicator: "Designing primers, please wait ... Time elapsed: 00 : 09".
- Progress steps: "Positioning LNA.", "Checking Tm.", "Checking secondary structure.", "Checking general specificity.".
- Text: "Please wait for the design." and "or".
- ⑦: Text input field for "Enter email to have a link to the result page sent by email:".
- "Send" button.

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Basket

Since you are located in a distributor country, Exiqon cannot offer you price information. To receive a quote, simply press the "Get quote" button below. You will be taken to a new page where you will be asked to submit your contact details. Your local distributor will then contact you with a quote.

test

Design ID: 520847-3
Product number: 309999
Description: ExiLERATE LNA qPCR custom primer set for 200 PCR reactions of 10µL
Note:

Remove

Quantity: 1

The following products are recommended with the items in your shopping basket:

Product no.: 303301
Product name: ExiLERATE LNA™ qPCR, cDNA Synthesis kit, 50 rxns
Description: ExiLERATE LNA™ qPCR, cDNA synthesis kit for reverse transcription of mRNA/lncRNA (50 rxns)

Please inquire

Product no.: 303402
Product name: ExiLERATE LNA™ qPCR, SYBR® Green master mix kit, 2.5ml
Description: ExiLERATE LNA™ qPCR, ExiLERATE SYBR® Green master mix, 2.5 ml (500 rxns of 10ul) and GAPDH endogenous control primer set

Please inquire

Product no.: 303410
Product name: ExiLERATE LNA™ qPCR, SYBR® Green master mix kit, 10ml
Description: ExiLERATE LNA™ qPCR, ExiLERATE SYBR® Green master mix, 10 ml (2000 rxns of 10ul) and GAPDH endogenous control primer set

Please inquire

Product no.: 309080
Product name: Random RT primer for LNA™ qPCR
Description: Random primer for use with Universal cDNA synthesis kit (product number 203301), 32 rxns

Please inquire

Press the 'Check out' button to continue the check out procedure online.

If you would rather like to email/fax your order, please follow these instructions:

- Press the 'Save quote - Order later' button.
- Email/Fax us this quote along with the following information:
 - Your name
 - Billing address
 - Shipping address
 - A copy of the purchase order
 - TAX/VAT ID number
- Email to: order@exiqon.com or use the following fax numbers:
 - North America: Fax: +1 781 376 4152
 - Outside North America: +45 45 66 18 88

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Contact us

In our effort to give you a rapid and fulfilling answer, we kindly ask you to fill in a few blanks below

Your inquiry is related to

- Technical Support
- Product Information
- Ordering
- Profiling Services
- Career in Exiqon
- Web & Marketing
- Sign up to Newsletter
- Other

- Sign up to our Newsletter and get research news, free downloads and offers 8-10 times a year. You can unsubscribe at anytime.

Please chose area of interest (click all that apply)

- microRNA Research
- mRNA Research
- Molecular Diagnostics
- Other

Please provide us with information about your inquiry. We will return to you as soon as the inquiry has been processed

Please send me a product quote.

Reference number: 45524

 [Click here to attach a file](#)

Reference number

Please state contact details

First name *

Last name *

E-mail *

Phone

Institution *

Address *

City *

Zip *

Country *

In which industry are you employed *

Are you currently engaged in microRNA research *

How did you first learn about Exiqon *

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