Recombinant Cas9 Protein GMP grade

Code No. T230 **Size:** **0.6 mg**
Conc.: **3 mg/ml**

Description:

Cas9 nuclease (*spCas9*) derived from *Streptococcus pyogenes* is widely used for gene editing in various cell types and organisms. CRISPR (Clustered Regularly Interspaced Short Palindromic Repeats)/Cas9 technology consists of two key components that form a complex: Cas9 nuclease and guide RNA (gRNA) that is a couple of small RNAs directing Cas9 nuclease to cleave a specific double-stranded genomic DNA in a sequence-specific manner^{1, 2, 3}.

Recombinant Cas9 Protein GMP grade is a recombinant wild-type *Streptococcus pyogenes* Cas9 nuclease containing nuclear-localization signal (NLS) at C-terminal expressed and purified from *E.coli*. The Cas9 protein solution has been verified to be well-tolerated by mammalian cells when electroporated or microinjected as a ribonucleoprotein complex (RNP) with a single guide RNA (sg RNA) for knockout experiments. This product is manufactured according to relevant GMP guidelines.

Storage : Store at or below -70°C

Molecular Mass:

159,665 (based on amino acid sequence, 1380 amino acids)

Source:

E. coli expressing recombinant wild-type *Streptococcus pyogenes* Cas9 nuclease with nuclear localization signal(NLS) of simian virus 40 at C-terminal.

Purity: ≥ 90% by HPLC

Packaging: 0.6 mg protein / 0.2 ml in a 1 ml cryovial

Application: gRNA-dependent double-stranded DNA cleavage

Precautions:

Avoid repeated freeze/thaw cycles.

* Takara Bio is not guaranteed the quality of the product after once opened by a customer. However, we have not found any obvious decline in performance after up to five(5) times freezing and thawing, if it is not opened.

Quality Statement:

1. This product is manufactured according to relevant GMP guidelines.
2. No animal- or human-derived components are used in the manufacture of this product.

References:

- 1) Jinek M, Chylinsky K, Fonfara I, Hauer M, Doudna J A, and Charpentier E. A programmable dual-RNA-guided DNA endonuclease in adaptive bacterial immunity. *Science*. (2012) **337 (6096)**: 816-821.
- 2) Liang X *et al.* Rapid and highly efficient mammalian cell engineering via Cas9 protein transfection. *J Biotechnol*. (2015) **208**: 44-53.
- 3) Sander J D and Joung J K. CRISPR-Cas9 systems for genomic editing, regulation and targeting. *Nat Biotechnol*. (2014) **32**: 347-355.

Related Products:

[Cas9 protein]
Guide-it™ Recombinant Cas9 (Electroporation-Ready)
(Cat. #632641/632640)

[Others]
Guide-it™ Cas9 Monoclonal Antibody (Clone TG8C1)
(Cat. #632628/632627)

Guide-it is a trademark of Takara Bio Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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● 起源
E. coli expressing recombinant wild-type *Streptococcus pyogenes* Cas9 nuclease with nuclear localization signal(NLS) of simian virus 40 at C-terminal.

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- 1) Jinek M, Chylinsky K, Fonfara I, Hauer M, Doudna J A, and Charpentier E. A programmable dual-RNA-guided DNA endonuclease in adaptive bacterial immunity. *Science*. (2012) **337 (6096)**: 816-821.
- 2) Liang X *et al.* Rapid and highly efficient mammalian cell engineering via Cas9 protein transfection. *J Biotechnol.* (2015) **208**: 44-53.
- 3) Sander J D and Joung J K. CRISPR-Cas9 systems for genomic editing, regulation and targeting. *Nat Biotechnol.* (2014) **32**: 347-355.

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[Cas9 タンパク質]
Guide-it™ Recombinant Cas9 (Electroporation-Ready)
(製品コード 632641/632640)

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v201910

Recombinant Cas9 Protein GMP grade

Experimental Example (Genome Editing of Mammalian Cells):

Reference: Guide-it Recombinant Cas9 (Electroporation-Ready) User Manual

[https://www.takarabio.com/assets/documents/User%20Manual/Guide-it%20Recombinant%20Cas9%20\(Electroporation-Ready\)%20User%20Manual_012317.pdf](https://www.takarabio.com/assets/documents/User%20Manual/Guide-it%20Recombinant%20Cas9%20(Electroporation-Ready)%20User%20Manual_012317.pdf)

Target cells / gene : 293T cell / CD81

RNP delivery method : Electroporation [Equipment: Neon Transfer System (Thermo Fisher Scientific)]

Preparation of RNP : 7.5 μ l solutions were prepared by mixed with 0.45 μ g of sgRNA, 0.75 μ g/0.25 μ l of this product and Resuspension Buffer R (or T), incubated at 37°C for 5 minutes and stand at 4°C.

RNP electroporation : Cells were harvested, washed with PBS and resuspended with Buffer R to be 2×10^7 cells /ml. RNPs were electroporated into cells using 7.5 μ l of RNP and 7.5 μ l of cells as prepared above according to equipment manual. Electroporation was conducted in the following three kinds of condition (N=3) (Table 1).

Table 1. Electroporation conditions

Condition	Pulse voltage (V)	Pulse interval (ms)	Pulse time
A	1,300	20.0	2
B	1,200	20.0	2
C	1,100	20.0	2

Efficiency of genome editing:

After the electroporation under the upper conditions, the cells were cultured for 7 days, and then reacted with PE-labeled Anti-CD81 antibody and measured the proportion of CD81 negative (knockout, KO) cells by FCM analysis.

Result:

Comparing Condition A to C, the highest Knockout (KO) efficiency of the CD81 gene (85.4%) was obtained at Condition A.

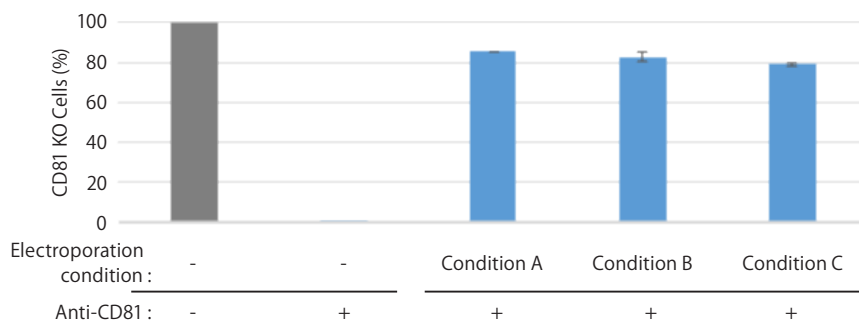
Figure 1. Knockout efficiency

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