# Brevibacillus Expression vector pNC-HisT DNA

Code No. HB121  
Size : 10  $\mu$ g (0.2 OD)

Shipping at  $-20^{\circ}\text{C}$   
Store at  $-20^{\circ}\text{C}$

\* 2 years from date of receipt under proper storage conditions.

Lot No.

Concentration : 0.2  $\mu$ g/ $\mu$ l

Volume : 50  $\mu$ l

Regarding protocol, please refer to the product manual supplied with this vector.

## Description :

pNC-HisT is a shuttle vector of *B. choshinensis* and *E. coli*, and this vector is a secretory expression vector. Expression plasmid can be constructed by using *E. coli* as the host, then transferred into *B. choshinensis* cells. The pNC-HisT contains P2 promoter derived from the host cell wall protein (HWP) as its expression promoter and does not work in *E. coli*. In *B. choshinensis* cells, P2 promoter works as very strong promoter, enabling efficient protein production.

His-Tag sequence is located at the downstream of the secretion signal peptide and Thrombin cleavage site is inserted so that the His-Tag can be removed from the fusion expression protein.

Form : 10 mM Tris-HCl, pH8.0  
1 mM EDTA

Base pairs : 5,260 bp


## Usage :

This is a vector for protein secretory expression using *Brevibacillus* Expression System (a shuttle vector between *B. choshinensis* and *E. coli*).

## Cloning sites of pNC-HisT DNA :

(Please refer to the other side of this sheet.)

## Map of pNC-HisT DNA :

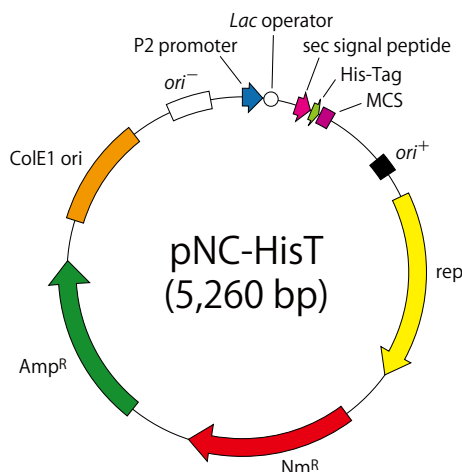## NOTICE TO PURCHASER: LIMITED LICENSE

### [L39] Brevibacillus Expression System

1. *Brevibacillus* Expression System-related products were developed and manufactured by Higeta Shoyu Co., Ltd. and sold by TAKARA BIO INC.
2. The use of this product is allowed for academic research only. You have to conclude a license agreement if you use this product or the products produced by using this product for commercial purpose. Contact us if you would like to ask any question about license etc. If any questions are received, either Higeta Shoyu Co., Ltd. or an agent for it will answer.  
TaKaRa Bio Inc.  
Tel +81-77-543-7247 Fax +81-77-543-9254
3. All rights of intellectual property related to this product are owned by Higeta Shoyu Co., Ltd. This product is covered by the following Japanese patents and corresponding international patent : Japanese patents, 3696322, 3734593, PCT/JP02/01038, PCT/JP2004/016912 (international application based on PCT treaty), 3433807.
4. This product, the components or the derivatives of this product, or the products produced by using these items can not be transferred (distributed freely or sold) to third party. (When you purchase this product, you submit the license agreement to us.)

## Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from TAKARA BIO INC. If you require licenses for other use, please contact us by phone at +81 77 543 7247 or from our website at [www.takara-bio.com](http://www.takara-bio.com). Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

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