pAAV-ZsGreen1 Vector

Code No. 6231

Size: 20 μ g

Conc.: 1.0 μ g/ μ l

* 2 years from date of receipt under proper storage conditions.

Regarding protocol, please refer to the product manual of AAVpro[®] Helper Free System (Cat. #6230/#6650/#6651/#6673)

Description:

pAAV-ZsGreen1 Vector is an adeno-associated virus (AAV) vector plasmid that has Inverted Terminal Repeat (ITR). This vector plasmid allows you to produce high titer recombinant AAV vector of serotype AAV1, AAV2, AAV5, or AAV6 together with AAVpro Helper Free System (AAV1) (Cat. #6673), AAVpro Helper Free System (AAV2) (Cat. #6230), AAVpro Helper Free System (AAV5) (Cat. #6650), or AAVpro Helper Free System (AAV6) (Cat. #6651).

This vector plasmid carries ZsGreen1 gene, a variant of the reef coral *Zoanthus sp.* green fluorescent protein, and allows ZsGreen1 to be used as an indicator of transfection efficiency and transduction efficiency.

Form:

10 mM Tris-HCl, pH 8.0
1 mM EDTA

Storage: -20°C

Preparation:


Purification of cccDNA by CsCl-EtBr ultracentrifugation and 0.22 μ m filtration

Length: 5,716 bp

Quality control:

1. Confirmed to maintain the 5' terminal sequence of ZsGreen1 by dideoxy sequencing method.
2. Shown to be cleaved at a single site with *Hinc* II and at four site with *Eco*O109 I.

Vector map of pAAV-ZsGreen1 Vector:

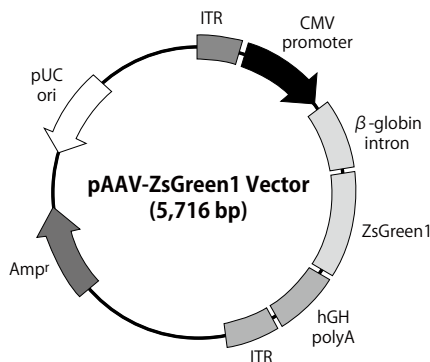Note:

- Please follow the guideline for experiments using recombinant DNA issued by the relevant authorities and the safety committee of your organization or your country in using this product.
- The adeno-associated virus (AAV) vector solution could contain potentially hazardous recombinant virus depending on an insert gene. Due caution must be exercised in this production and handling of recombinant AAV vector.
- Takara Bio is not liable for any accident or damage caused by the use of this product.

AAVpro is a registered trademark of Takara Bio Inc.

Living Colors is a registered trademark of Takara Bio USA, Inc.

Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6973 or from our website at www.takara-bio.com.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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