# pAAV8-ZsGreen1 One-Vector

**Code No. 6268**

**Size: 500  $\mu$ l x 2**

**Conc.: 1.0  $\mu$ g/ $\mu$ l**

\* 2 years from date of receipt under proper storage conditions.

## Description:

pAAV8-ZsGreen1 One-vector is a single plasmid that incorporates all three elements required for AAV vector production: the transfer plasmid (pAAV-ZsGreen1), the Rep/Cap plasmid (pRC8), and the adenoviral helper plasmid (pHelper).

The ZsGreen1 reporter gene is located between the ITRs and is driven by the CMV promoter, allowing efficient expression of ZsGreen1 in target cells following AAV transduction.

This plasmid enables the production of AAV particles expressing ZsGreen1 without the need for co-transfection of multiple plasmids.

The ZsGreen1 gene can be removed by *Not* I digestion, allowing users to replace it with their gene of interest. For cloning, it strongly recommends using In-Fusion® Snap Assembly Master Mix (Cat. #638943/638944/638947 to 638949).

**Note:** Due to the packaging size limitation of AAV particles, the inserted gene should be no larger than 2.8 kb.

**Form:** 10 mM Tris-HCl, pH 8.0  
1 mM EDTA

**Storage:** -20°C

## Preparation:


Purification of cccDNA by CsCl-EtBr ultracentrifugation and 0.22  $\mu$ m filtration.

**Length:** 19,282 bp

## Quality Control Data:

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

## Vector Map of pAAV8-ZsGreen1 One-Vector:

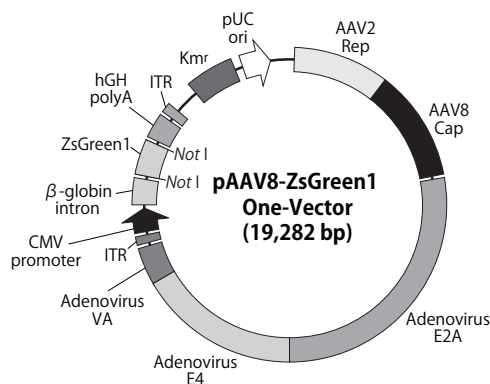## Precautions for Use:

- Please follow the guideline for experiments using recombinant DNA issued by the relevant authorities and the safety committee of your organization or your country in using this product.
- The adeno-associated virus (AAV) vector solution could contain potentially hazardous recombinant virus depending on an insert gene. Due caution must be exercised in this production and handling of recombinant AAV vector.
- Takara Bio is not liable for any accident or damage caused by the use of this product.

In-Fusion is a registered trademark of Takara Bio USA, Inc.

## Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us from our website at [www.takarabio.com](http://www.takarabio.com).

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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