SeV Primer/Probe Mix with Standard

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Cat. No. QS-0100-200

I. Product Description

The SeV Primer/Probe Mix enables highly sensitive detection of Sendai virus (SeV), making it ideal for confirming viral presence in infected cells. Additionally, SeV genome can be quantified using the Standard and Standard Dilution Buffer included in the kit.

The assay can be performed as either a one-step assay or two-step procedure. Cell lysate or extracted total RNA can be used as templates.

Preparation example for one reaction

One Step PrimeScript™ III RT-qPCR Mix (2×)	10μL
SeV Primer/Probe Mix (×10)	2μL
RNase Free Water	6μL
Template	(2μL)
<hr/>	
Total	20μL

II. Product Contents

Product Component	Quantity	Cat No.
SeV Primer/Probe Mix (×10)	400μl	Q-0120
SeV Standard (2.0E+06 copies/μl)	120μl	QD-0103
Standard Dilution Buffer (×200)	50μl	QB-0105

(Probe: 5'FAM-3'MGB-Eclipse label)

III. Real-time PCR

1. Preparation of Standard Dilution Buffer

Prepare a 1×Standard Dilution Buffer by diluting the 200 × Standard Dilution Buffer 200-fold with RNase-Free Water.

2. Preparation of SeV Standard

Prepare 10-fold serial dilutions of SeV Standard using the SeV Standard stock (2×10⁶ copies/μL) and Standard Dilution Buffer as follows:

(Use 2μL of template for each real-time PCR reaction)

- (1) 2×10⁶ copies/μL
[SeV Standard stock]
- (2) 2×10⁵ copies/μL
[2μL SeV Standard stock+18μL Standard Dilution Buffer]
- (3) 2×10⁴ copies/μL
[2μL of dilution (2)+18μL Standard Dilution Buffer]
- (4) 2×10³ copies/μL
[2μL of dilution (3)+18μL Standard Dilution Buffer]
- (5) 2×10² copies/μL
[2μL of dilution (4)+18μL Standard Dilution Buffer]
- (6) 2×10¹ copies/μL
[2μL of dilution (5)+18μL Standard Dilution Buffer]

3. PCR Setup

Perform real-time PCR according to the instructions provided with the One Step PrimeScript™ III RT-qPCR Mix.

4. qPCR Cycling Conditions

The following conditions are recommended for the LightCycler® 480 System II. When using different real-time PCR instruments, optimization of cycling conditions may be required.

Step	Temperature	Duration	Cycle(s)
Reverse Transcription	52°C	5 min	1
	95°C	10 sec	
2 Step PCR	95°C	5 sec	55
	60°C	30 sec	

Fluorescence detection: FAM

IV. Example of Results

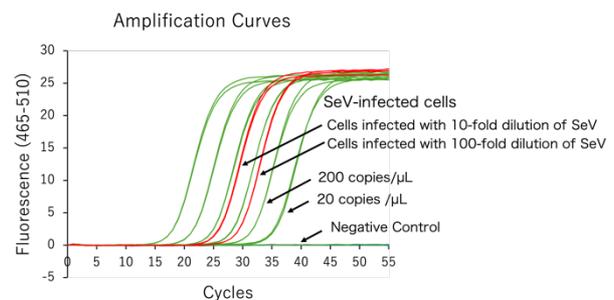
Figure 1: Example of SeV real-time PCR using SeV Primer/Probe Mix with Standard

V. Storage

All product components should be stored at -20°C and used within 6 months after delivery.

VI. Product use restriction

· This product is for research use only and should not be used for diagnostic or therapeutic purposes.