BspQ I GMP grade

G C T C T T C N |
C G A G A A G N N N N

Code No. T2451S001	Size: 20 KU
	Conc.: 20 U/ μ l
Code No. T2451S010	Size: 200 KU
	Conc.: 20 U/ μ l
Code No. T2451S100	Size: 2,000 KU
	Conc.: 20 U/ μ l

Description:

This enzyme has an equivalent performance to BspQ I, HQ (Cat. #1228A), which recognizes asymmetric specific double-stranded DNA sequences and cleaves both strands outside the recognition sequence (Type IIS restriction enzyme). Therefore, in the preparation of linearized plasmids for *in vitro* transcription (IVT), introducing the recognition site of this enzyme immediately after the Poly(A) sequence of the target mRNA makes it possible to design mRNAs that do not have an extra sequence after the Poly(A) sequence.

Quality Statement:

1. This product is manufactured in accordance with relevant GMP guidelines. For more information, please refer to our website.
2. No animal- or human-derived components are used in the manufacture of this product.
3. This product uses primary materials that do not contain β -lactam compounds.

Quality Control Data:

Please see the Certificate of Analysis (CoA) for each lot. You can download the CoA on Takara Bio website.

Storage Buffer:

20 mM	Tris-HCl, pH 7.5
500 mM	KCl
0.1 mM	EDTA
1 mM	DTT
0.1%	Tween 20
50%	Glycerol

Storage: $-20 \pm 5^\circ\text{C}$

Source:

Escherichia coli carrying the plasmid containing the gene for BspQ I

Unit Definition:

One unit is defined as the amount of enzyme that cleaves 10 pmol of fluorescence-labeled probes in 1 hour at 50°C in a 50 μ l reaction.

Application:

Linearization of template plasmid used for mRNA synthesis

Precautions for Use:

Do not mix the enzyme vigorously.

General Reaction Mixture and Incubation Time:

BspQ I GMP grade*1	1 μ l
10X BspQ I Buffer GMP grade*2	10 μ l
Substrate DNA	$\leq 10 \mu\text{g}$
Nuclease-free water	up to 100 μ l

↓

Incubate at 50°C for 1 hour.

*1 Add this to the reaction last.

*2 Cat. #T2461S010/S100 (sold separately)

Reaction Temperature: 50°C

Note: The enzyme activity at 37°C decreases to about 30% of its activity at 50°C.

Heat Inactivation: 70°C for 20 min

Effect of DNA Methylation:

This enzyme is not sensitive to Dam, Dcm, and CpG methylation.

Star Activity:

Star activity may result from an excess incubation (e.g., > 4 hours).

Related Products:

10X BspQ I Buffer GMP grade (Cat. #T2461S010/S100)

T7 RNA Polymerase GMP grade (Cat. #T2411S001/S010/S100)

PrimeCap® T7 RNA Polymerase (low dsRNA) GMP grade

(Cat. #T2422S001/S010/S100)

10X IVT Reaction Buffer GMP grade (Cat. #T2401S010/S100/S500)

Recombinant RNase Inhibitor GMP grade (Cat. #T2431S001/S010/S100)

PrimeCap is a registered trademark of Takara Bio Inc.

Note

This product is not intended for humans or animals *in vivo* applications. Our products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without our prior written approval.

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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