# EDTA Buffer Powder, pH 8.0

**Code No. T9191**      **Size: 5 pouches**  
**(for 2.5 L)**

## Description:

EDTA (ethylene-diamine-tetraacetic acid) is a chelating agent to sequester divalent metal ions such as calcium and magnesium. This ability prevents DNA and RNA degradation as metal-dependent enzymes acting as nucleases become deactivated.

This product is a powder for preparing EDTA solution which can easily be prepared by dissolving it with H<sub>2</sub>O. One pouch is for preparing 500 ml of 0.5 M EDTA (pH 8.0).

## Storage:

Room temperature at desiccated condition.

## Specifications:

Format: Powder  
Concentration: 0.5 M EDTA  
Volume: For 500 ml  
pH: 8.0 ± 0.05 at 25°C

## Preparation for use:

Transfer one pouch of EDTA Buffer Powder, pH 8.0 in a flask or beaker. Add 300 ml of distilled water or deionized water and stir the solution for a few minutes. Adjust the volume to 500 ml, and stir until complete dissolution.

## Applications:

- Inhibitor and stop solution for metal-dependent enzymatic reaction
- Preparation of TE Buffer

### Note

This product is for research use only. It is not intended for use in therapeutic or diagnostic procedures for humans or animals. Also, do not use this product as food, cosmetic, or household item, etc. Takara products may not be resold or transferred, modified for resale or transfer, or used to manufacture commercial products without written approval from Takara Bio Inc.

If you require licenses for other use, please contact us by phone at +81 77 565 6972 or from our website at [www.takarabio.com](http://www.takarabio.com).

Your use of this product is also subject to compliance with any applicable licensing requirements described on the product web page. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.

All trademarks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions.

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