

# Fastest, Easiest Adenoviral System Ever

## Adeno-X™ Adenoviral System 3

- Cloning into adenovirus is as simple and rapid as into any plasmid
- Clone directly into the pAdenoX vector, no shuttle vector required
- Transduce dividing and non-dividing cells
- Very high titer, high expression, broad host range
- Highly flexible formats


The **Adeno-X Adenoviral System 3** is the most advanced commercially available adenoviral gene delivery system—providing by far the simplest, fastest, and most efficient method for constructing recombinant adenoviral vectors.

### Why choose adenoviral gene delivery?

Adenoviral gene transfer is one of the most reliable methods for introducing genes into mammalian cells. Because infection by adenovirus is not cell-cycle dependent, you can deliver your gene to primary as well as transformed cell lines. Adenoviruses are ideal tools for protein production in mammalian cells because following infection, your target gene is transiently expressed at very high levels since many cells receive multiple copies of the recombinant genome. Additionally, adenoviruses are capable of infecting a wide variety of proliferating and quiescent cell types from many different animal species including humans, non-human primates, pigs, rodents, mice, and rabbits.

### Clone into adenovirus just like any other plasmid

Until now the main drawback of commercially supplied adenoviral vector systems has been the need to use complex cloning procedures to overcome the difficulties with cloning into large (~34 kb) plasmids. Procedures have included precloning into shuttle vectors and subcloning through multiple steps and multiple different strains of *E. coli*, all of which increase hands-on time and gives more room for error. At Clontech, our Adeno-X virologists thought “wouldn't it be great if you could clone directly into the adenoviral plasmid just like any plasmid?” They then harnessed the power of In-Fusion® HD cloning technology to make this happen.

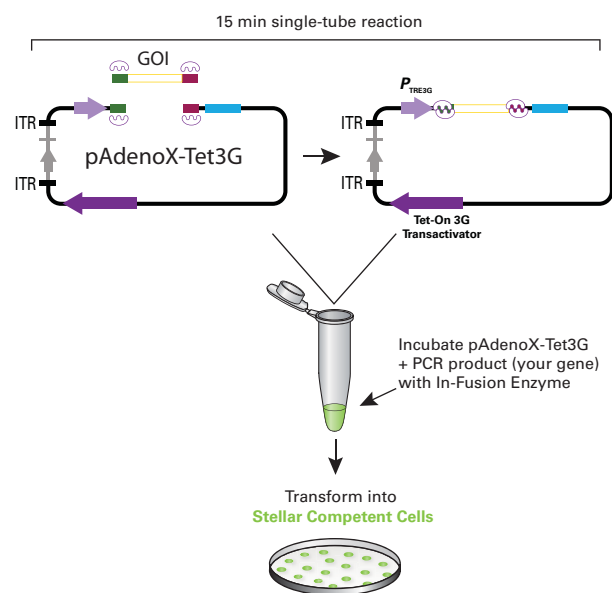

### Adeno-X System 3 Features

Feature	Description
No shuttle vector required	Clone directly into the adenoviral vector using In-Fusion HD
Easy to use	As easy as any plasmid cloning system
Fast	2-3 days for cloning (others take at least 8 days)
Super high-efficiency cloning	100s of colonies, 9/10 clones are correct
Highly flexible formats	Use an existing expression cassette or create one without additional subcloning
Best technologies offered	Tet-On® 3G Inducible Expression, fluorescent reporters, multiple fragment cloning

**Clontech Laboratories, Inc.**  
A Takara Bio Company  
[www.clontech.com](http://www.clontech.com)


United States/Canada: +1.800.662.2566 • Asia Pacific: +1.650.919.7300 • Europe: +33.(0)1.3904.6880 • Japan: +81.(0)77.543.6116  
For Research Use Only. Not for use in diagnostic or therapeutic procedures. Not for resale. Clontech, the Clontech logo, Adeno-X, In-Fusion, Stellar and Tet-On are trademarks of Clontech Laboratories, Inc. All other marks are the property of their respective owners. Certain trademarks may not be registered in all jurisdictions. ©2011 Clontech Laboratories, Inc. AN143880 US (633288)



# Fastest, Easiest Adenoviral System Ever ...continued

## There is no simpler adenoviral expression system

An overview of the procedure for creating recombinant adenovirus using the Adeno-X Adenoviral System 3 is shown in **Figure 1**. The system relies upon the ability of the In-Fusion HD enzyme to precisely recognize and fuse 15 bp of homology between two linear DNA molecules. To generate short regions of homology to the prelinearized pAdeno-X vector, simply add 15 bp of additional sequence to the primers you use to PCR amplify your gene of interest. Then, combine the DNA together with the In-Fusion HD enzyme in a 15 min reaction and transform Stellar Competent Cells. Cloning is always directional and 90% of the clones contain the correct insert. The kit includes a control cloning fragment, primers for colony PCR verification of positive clones and a NucleoBond Xtra kit for transfection-grade plasmid purification.

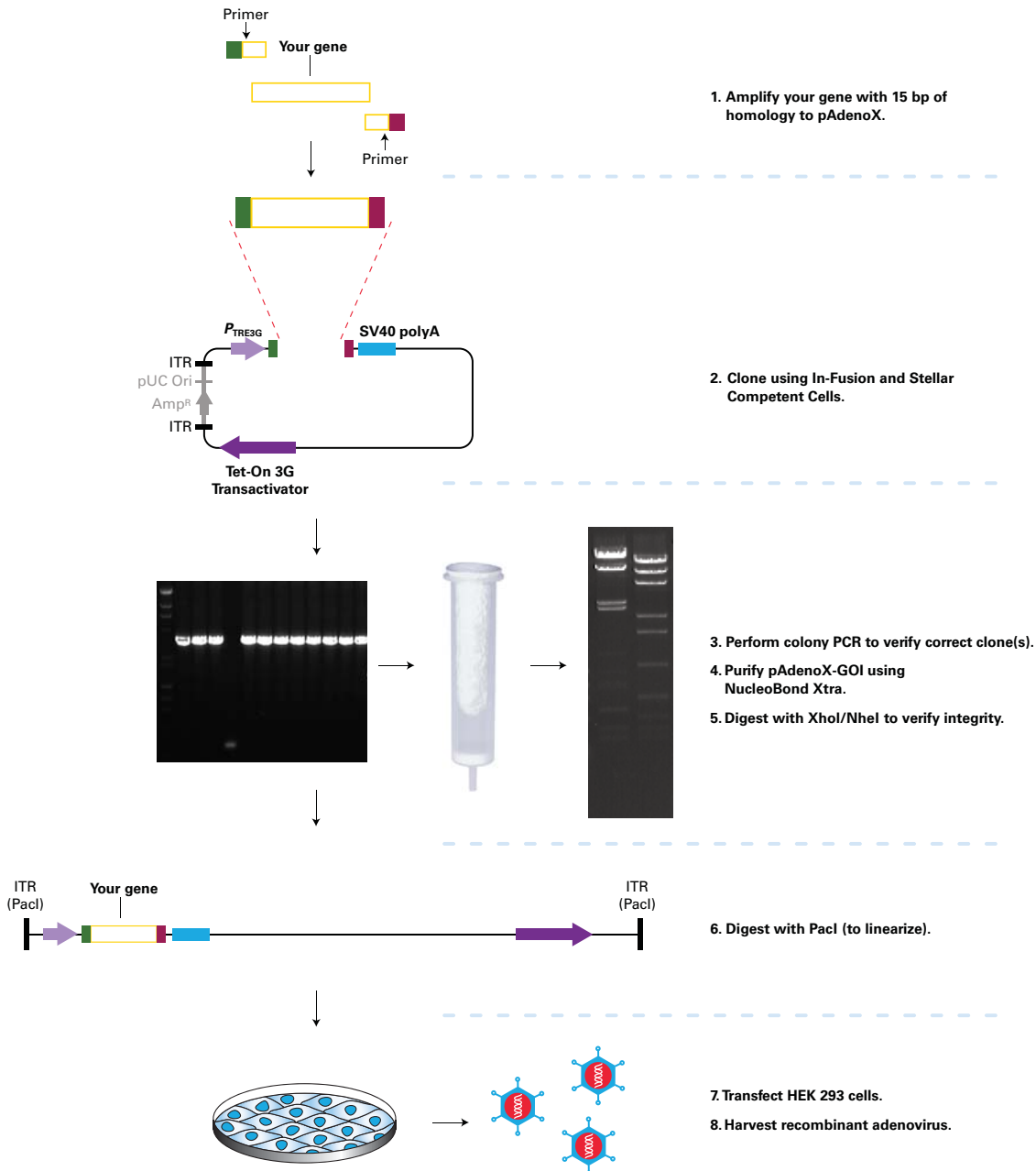

**Figure 1. Constructing recombinant adenovirus with In-Fusion technology.** DNA sequences can be rapidly transferred as PCR products to any pAdenoX vector using the In-Fusion cloning method. In this example, your gene of interest is amplified with 15 bp extensions that are homologous to the ends of the linearized adenoviral vector. The PCR product is then purified and mixed with the linearized adenoviral vector of choice in the In-Fusion reaction. Following the reaction, a portion of the mixture is transformed into Stellar™ Competent Cells and screened. Once a PCR-positive clone is identified, the recombinant pAdenoX vector is amplified, purified, and subsequently linearized with the restriction enzyme PacI, then transfected into HEK 293 cells for viral rescue and amplification.

# Fastest, Easiest Adenoviral System Ever ...continued

## Multiple formats are available

Adeno-X Adenoviral System 3 is available in seven formats, including the most advanced tetracycline inducible expression system, constitutive expression systems with or without fluorescent reporters, and universal systems that allow you to clone and express any entire expression cassette of your choice (**Table I**).

Table I: Adeno-X Adenoviral System 3 Formats			
Cat. No.	Product	Description	Vector Map
631180	Adeno-X Adenoviral System 3 (Tet-On 3G Inducible)	<ul style="list-style-type: none"> <li>Tightly-controlled, doxycycline-inducible expression system</li> </ul>	
632269	Adeno-X Adenoviral System 3 (CMV)	<ul style="list-style-type: none"> <li>Constitutive expression from a CMV promoter</li> </ul>	
632268	Adeno-X Adenoviral System 3 (CMV, Red)	<ul style="list-style-type: none"> <li>Constitutive expression from a CMV promoter</li> <li>Red fluorescent protein to easily monitor transfection and transduction</li> </ul>	
632267	Adeno-X Adenoviral System 3 (CMV, Green)	<ul style="list-style-type: none"> <li>Constitutive expression from a CMV promoter</li> <li>Green fluorescent protein to easily monitor transfection and transduction</li> </ul>	
632266	Adeno-X Adenoviral System 3 (Universal)	<ul style="list-style-type: none"> <li>Use any promoter and any polyA sequence</li> <li>Ideal for tissue-specific expression or expression of shRNA or miRNA</li> </ul>	
632265	Adeno-X Adenoviral System 3 (Universal, Red)	<ul style="list-style-type: none"> <li>Use any promoter and any polyA sequence</li> <li>Ideal for tissue-specific expression or expression of shRNA or miRNA</li> <li>Red fluorescent protein to easily monitor transfection and transduction</li> </ul>	
632264	Adeno-X Adenoviral System 3 (Universal, Green)	<ul style="list-style-type: none"> <li>Use any promoter and any polyA sequence</li> <li>Ideal for tissue-specific expression or expression of shRNA or miRNA</li> <li>Green fluorescent protein to easily monitor transfection and transduction</li> </ul>	


# Fastest, Easiest Adenoviral System Ever...continued

## Tetracycline inducible expression

When you clone your gene into pAdenoX-Tet3G you are creating a system with the tightest and most sensitive control of gene expression. Tightly-controlled, doxycycline-induced expression is as easy as constitutive expression since the Tet-On 3G transactivator protein and the  $P_{TRE3G}$  controlled gene of interest are contained on the same adenoviral vector. Up to 3000-fold induction can be achieved using this system (Figure 3).


## Clone any expression cassette into the universal vectors

You are not limited to using a CMV expression system—we have created universal systems with vectors that lack a promoter and polyA signal in the cloning site. Simply amplify an entire expression cassette (from promoter to polyA) from a pre-existing construct and clone using In-Fusion HD (Figure 4, Panel A). Universal systems can be used for expression from alternative promoters that are more suitable to your target cell type such as EF1 alpha or tissue-specific promoters. Alternatively, you may wish to transfer your shRNA or miRNA expression cassette from a pre-existing plasmid to one of the universal pAdenoX vectors in order to create a high efficiency RNAi delivery system. Even if your expression cassette does not exist you can create one using multiple fragment cloning (Figure 4, Panel B), with only a small loss in cloning efficiency (Table II).

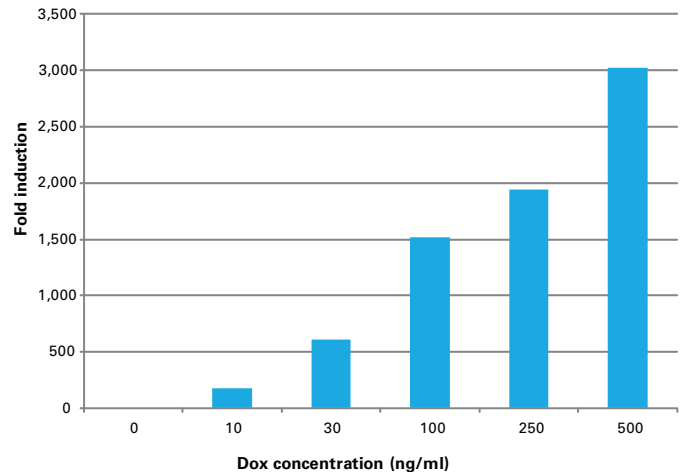


**Figure 3. The Adeno-X Tet-On 3G Systems generate very high-fold induction, with up to 3000-fold difference between induced and uninduced states.** Using equal amounts of high-titer supernatants, HeLa cells cultured at the indicated concentrations of Dox were infected with Adeno-X Tet-On 3G Luciferase virus. Cultures were harvested and assayed for luciferase activity.

### A Complete Expression cassette cloned into a Universal System

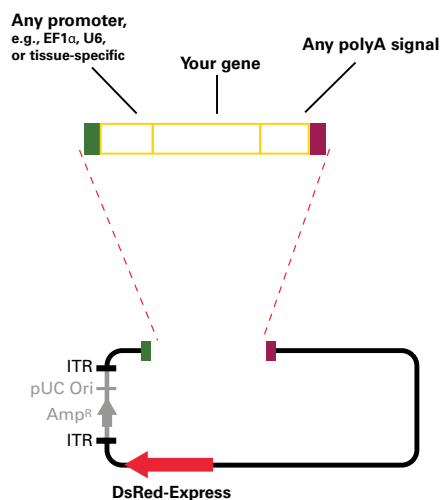

### B Multiple fragment cloning in the Universal Red System

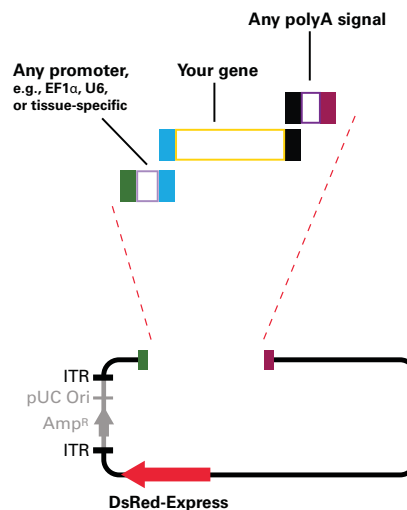

**Figure 4. The Universal Adeno-X Expression Systems contain vectors that lack a promoter and polyA signal in the cloning site.** You can either clone an expression cassette from a preexisting construct into the vector (Panel A) or create a new one using multiple fragment In-Fusion HD cloning (Panel B).

**Table II: Multiple Fragment Cloning Efficiency with the Adeno-X System 3**

	Number of colonies (1/10th plated)	Percentage correct
pAdenoX + 1 insert	200-300	90%
pAdenoX + 2 inserts	200-300	60%
pAdenoX + 3 inserts	25-40	25%

# Fastest, Easiest Adenoviral System Ever ...continued

## Very high titer, easily amplified, very stable

Recombinant adenoviruses such as Adeno-X are lytic only in packaging cells that provide the essential E1 protein *in trans* (such as HEK 293 cells). This lytic mechanism of amplification means that virus particles produced by one cell can reinfect adjacent packaging cells to produce a cascade of virus production and ultimately far higher titers ( $>10^9$  IFU/ml) than can be achieved with any recombinant lentivirus system (**Table III**). Moreover, it is very simple to reamplify and make more virus particles. Unlike lentivirus production, there is no need to optimize transfection conditions, you simply reinfect HEK 293 cells with your existing adenovirus prep and wait for the packaging cells to produce more virus. Adenovirus can be stored in high titer aliquots frozen for long-term studies.

**Table III : Adenoviral Gene Delivery vs Lentiviral Gene Delivery**

	Lentivirus	Adenovirus
Infects many different human cell types	Yes	Yes
Infects both dividing and non-dividing cells	Yes	Yes
Non-integrating virus	No	Yes
High level of protein expression (up to 10-20% total protein)	No	Yes
Ability to accommodate long inserts (up to 8 kb)	No	Yes
Easy to scale-up/amplify	No	Yes
Easy to get titers $>10^9$ IFU/ml	No	Yes
Easy to get a multiplicity of infection $>25$ copies per cell	No	Yes

Product	Size	Cat. No.
Adeno-X Adenoviral System 3 (Tet-On 3G Inducible)	each	631180
Adeno-X Adenoviral System 3 (CMV)	each	632269
Adeno-X Adenoviral System 3 (CMV, Red)	each	632268
Adeno-X Adenoviral System 3 (CMV, Green)	each	632267
Adeno-X Adenoviral System 3 (Universal)	each	632266
Adeno-X Adenoviral System 3 (Universal, Red)	each	632265
Adeno-X Adenoviral System 3 (Universal, Green)	each	632264

### Notice to Purchaser

Your use of these products and technologies is subject to compliance with any applicable licensing requirements described on the product's web page at <http://www.clontech.com>. It is your responsibility to review, understand and adhere to any restrictions imposed by such statements.